



TRANSCCELL TECHNOLOGY, INC.

PBS

Precision Balance

Operation & Calibration Manual

Revision 2.3
MAY 19, 2003

© 2002 Transcell Technology, Inc.

Contents subject to change without notice.

Transcell Technology, Inc.
975 Deerfield Parkway
Buffalo Grove, IL 60089
Tel (847) 419-9180
Fax (847) 419-1515
E-mail: transcell@transcell.net
Web: www.transcell.net

Electromagnetic Compatibility Statement for Europe

PRODUCT	PBS Series	Directives with which this equipment complies:
Harmonized Standards applied:		EMC 89/336/EEC EMC Directive
		EMC 92/31/EEC EMC Directive
		EMC 93/68/EEC EMC Directive
EN 55022:1998 (Class B)		
EN 61000-3-2:1995 + A1:1998 + A2:1998		
EN 61000-3-3:1995		
EN 55024:1998		EN
61000-4-2:1995 + A1:1998		
EN 61000-4-3:1996		
EN 61000-4-4:1995		
EN 61000-4-5:1995		
EN 61000-4-6:1996		
EN 61000-4-8:1993		
EN 61000-4-11:1994		
Year in which CE mark affixed	2001	Date of issue October 2001

Electromagnetic Compatibility Statement for North America

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his/her own expense.

TABLE OF CONTENTS

	<u>Page</u>
Chapter 1: Introduction to the Transcell PBS Series Precision Balance	5
Chapter 2: Getting Started.....	6
Chapter 3: Operation	7
3.1 Display	7
3.1.1 Liquid Crystal Display (LCD)	7
3.2 Keyboard.....	8
3.2.1 Function Keys	8
3.3 General Scale Operation	8
3.3.1 Weighing an Item.....	8
3.3.2 Taring an Item.....	9
3.3.3 Clearing a Tare	9
Chapter 4: Advanced Features and Operation.....	10
4.1 Piece Counting.....	10
4.2 Clearing the Piece Count	11
4.3 Memory Accumulator	11
4.3.1 Displaying the Memory Accumulator	11
4.3.2 Adding to the Memory Accumulator.....	11
4.3.3 Clearing the Memory Accumulator.....	11
4.4 Percentage Mode.....	11
4.5 Display APW Mode.....	12
Chapter 5: Scale Configuration.....	13
5.1 Scale Configuration Overview.....	13
5.2 User (“A”) Menu	13
5.2.1 Entering the User Menu	13
5.2.2 Navigating in the User Menu.....	13
5.2.3 Notes on the User Menu	14
5.2.4 Units Selection Procedure (A6).....	15
5.2.5 Custom Unit Procedure (A7).....	16
5.2.6 Exiting the User Menu.....	16
5.3 Factory Setup (“F”) Menu.....	16
6.1 Scale Calibration Procedure	19
6.2 Calibration adjustment	20

Appendix A: Specifications	21
Appendix B: Serial Port Information	23
B.1 Serial Port Hardware	23
B.2 Serial Port Modes	23
B.2.1 Demand Mode	23
B.2.2 Continuous Mode	25
B.2.3 Recognized Host Commands	26
B.3 Transmit APW Mode	26
Appendix C: Error Messages	27
C.1 Error Messages	27
C.1.1 Operator Errors	27

LIST OF FIGURES

1-1 PBS Series Front Panel	5
2-1 PBS Back Panel	6
3-1 PBS Display Detail	7
3-2 Function Keys Layout	8
5-1 User Menu Key Assignments	13
5-2 User Menu Chart	14
6-1 Calibration Mode Key Assignments	19
B-1 Pin Assignments for DSUB9 serial port connector	22
B-2 Cable Diagram for Scale to PC	22
B-3 Cable Diagram for Scale to Printer	22
B-4 Demand Mode Output Format for Print Command	23
B-5 Demand Mode Output Format for M+ Command	23
B-6 Demand Mode Output Format for MR Command	24
B-7 Continuous Mode Output Format	24

LIST OF TABLES

1-1 PBS Series Product Matrix	5
3-1 PBS Series Annunciator Definitions	7
4-1 PBS Piece Counting Sampling Limits	10
4-3 PBS Percentage Sampling Limits	12
5-1 User Menu Chart	14
6-1 PBS Calibration Test Weight	19

CHAPTER 1: INTRODUCTION TO THE TRANSCELL PBS SERIES PRECISION BALANCE

The PBS Series Precision Balance, available in several capacities, is a large capacity, affordable balance designed for the classroom, laboratory or industrial applications. The PBS features 1:30,000 displayed resolution, a large, high contrast LCD display and digital calibration. Weighing modes include gram (kilogram), carat, avoirdupois pound and troy ounce. In addition, you may program one additional custom unit. A “Plus” model is available that features 1:50,000 displayed resolution.

The scale is available in three weight capacities as illustrated in Table 1-1. The “Plus” model matrix is shown in Table 1-2.

Prior to using the scale, please read this user’s guide carefully and completely. Store the manual in a safe and convenient place so it will be available if you have questions concerning the operation of the scale.

MODEL	CAPACITY / GRADUATION
PBS-300	300 x 0.01 g
PBS-600	600 x 0.02 g
PBS-1500	1500 x 0.05 g
PBS-3000	3000 x 0.1 g

Table 1-1: PBS Series Product Matrix

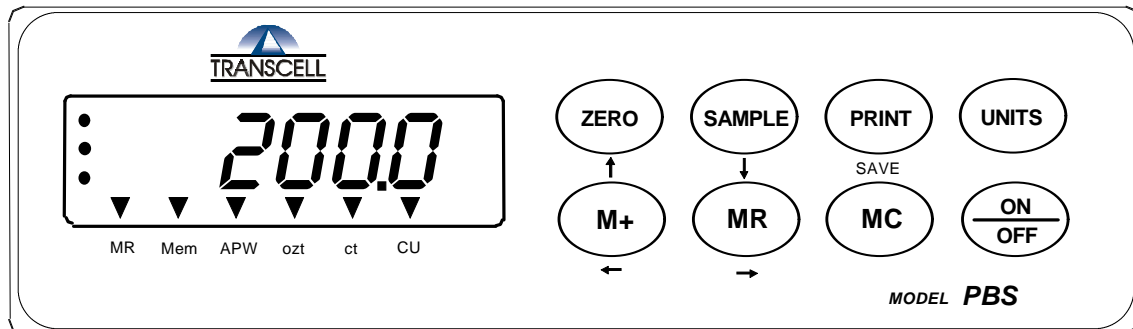


Figure 1-1: PBS Series Front Panel

CHAPTER 2: GETTING STARTED

After unpacking the scale, a small amount of preparation is required before the scale can be used. Please refer to Figure 2-1 below as needed.

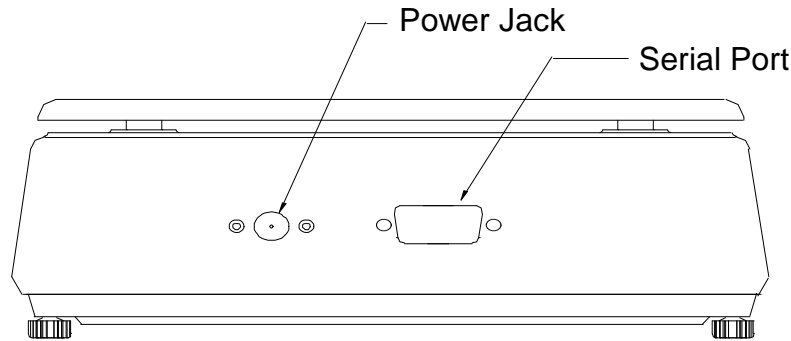


Figure 2-1: PBS Back Panel

- Step 1. Position the scale in its area of intended use. Observe the following guidelines for suitable location.
1. Choose a firm, stable floor or table.
 2. Do not share an AC outlet with electrical noise producing equipment, such as refrigeration units. This includes products with electrical motors and/or relays.
 3. Do not place the scale in an area with changing ambient temperature and/or high humidity.
 4. Do not place the scale in an area prone to exposure to direct sunlight, wind, or dust.
 5. Do not place the scale in an area with vibrating equipment.
- Step 2. Install the AC Adapter.
1. After placing the scale in its area of use, locate the AC Adapter.
 2. Connect the female end of the AC Adapter to the connector on the rear of scale, and then plug the adapter into an AC outlet. ***Make sure that the AC voltage appearing at the wall outlet matches the input voltage marked on the AC adapter.***
- NOTE:** Six AA size alkaline batteries may also be used to power the scale. The battery holder is located underneath the scale. Please observe the indicated polarity when installing the batteries.
- Step 3. If applicable, install the serial printer.
1. Connect the printer to the PBS's interface port using the optional serial cable.
 2. Configure the printer's communication parameters as detailed in Chapter 5.
- Step 4. Push the ON/OFF key on the front panel to begin use.

CHAPTER 3: OPERATION

3.1 DISPLAY

The Model PBS series scale utilizes a 6 character LCD (Liquid Crystal Display) to display the weight and system information.

3.1.1 LIQUID CRYSTAL DISPLAY (LCD)

Figure 3-1 shows the display detail of the PBS Series. The display consists of 6 numeric characters, weight units and system status flags. Table 3-1 lists the various annunciators you may see and their meanings.

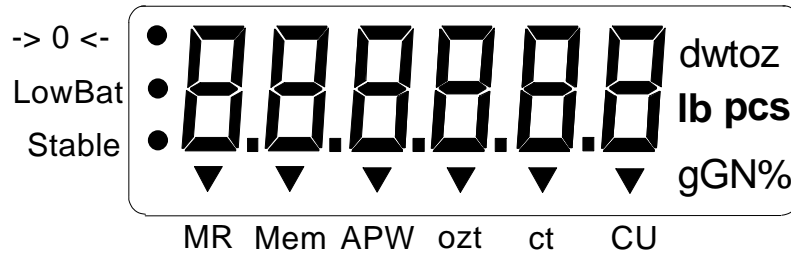


FIGURE 3-1: PBS Display Detail

Annunciator	MEANING
-> 0 <-	Center of Zero annunciator. Indicates that the scale is at gross zero.
Low Bat	Indicates a low battery condition. Check / replace the batteries.
Stable	This light is on whenever the scale is stable.
MR	Indicates that the scale is briefly displaying the contents of the memory accumulator.
Mem	Indicates that there is a value greater than zero in the memory accumulator.
APW	Indicates that the scale is displaying the APW (Average Piece Weight) in grams of the items you are counting.
ozt	Indicates that the scale is currently weighing in troy ounces.
ct	Indicates that the scale is currently weighing in carats.
CU	Indicates that the scale is currently weighing in your programmed custom unit.
dwt, oz, lb, g, GN	Indicates the unit of measure the scale is currently weighing in.
pcs, %	Indicates that the scale is currently in piece count (pcs) or percentage mode (%).

TABLE 3-1: PBS Series Annunciator Definitions

3.2 KEYBOARD

The keyboard is composed of eight function keys. Refer to Figure 3-2 for the overall layout and key locations.

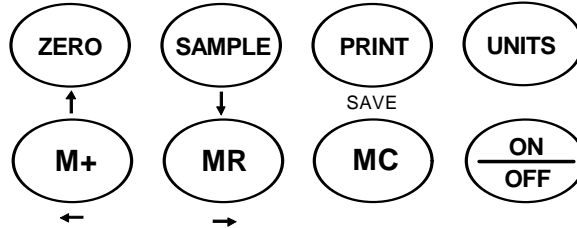


FIGURE 3-2: Function Keys Layout

3.2.1 FUNCTION KEYS

Zero - This key sets the scale to display zero.

Sample – This key puts the scale into sampling mode, which is used for piece counting. See Section 4.1 for more information.

Print - This key is used to send weight information out to the serial interface port. The scale cannot be in motion or in an overload condition.

Units – This key toggles the scale among the available weight units. The available weight units are lb, oz, g, pieces, %, GN, dwt, ct, ozt and custom unit. To see how to turn some units off, see Section 5.2.4.

M+ – Pressing this key adds the displayed weight or pieces to the memory accumulator. See Section 4.3 for more information.

MR – Pressing this key briefly displays the contents of the memory accumulator but only if the Mem annunciator is lit. See Section 4.3 for more information

MC – Pressing this key clears the memory accumulator. See Section 4.3 for more information.

On/Off - This key toggles the scale ON and OFF.

3.3 GENERAL SCALE OPERATION

3.3.1 WEIGHING AN ITEM

1. Select the desired weighing unit by pressing the UNITS key until that unit is indicated on the display.
2. If necessary, press the ZERO key to obtain a weight reading of zero.
3. Place the object to be weighed on the scale's platter and allow the weight indication to stabilize. If the item weight exceeds the scale's weight capacity, it displays "ooooo".
4. Read the weight shown on the display.

3.3.2 TARING AN ITEM

To weigh an item in a container, the weight of that container must first be subtracted from the overall weight to obtain an accurate weight reading. This is known as taring.

1. Select the desired weighing unit by pressing the UNITS key until that unit is indicated on the display.
2. Place the empty container on the scale's platter and allow the weight indication to stabilize.
3. Press the ZERO key.
4. Place the material to be weighed in the container and allow the weight indication to stabilize.
5. Read the weight shown on the display.

3.3.3 CLEARING A TARE

1. To clear a tare, take everything off of the scale's platter and press the ZERO key.

CHAPTER 4: ADVANCED FEATURES AND OPERATION

4.1 PIECE COUNTING

This mode is used to indicate the number of pieces of an item you have placed on the scale's platform and is accessed by pressing the SAMPLE key. To ensure accuracy, the parts you are counting must be consistent in weight.

The scale uses the sampling method to determine the average piece weight (APW) of the items you wish to count. When sampling items, always count the parts in your hand and place them on the platform all at once. If the APW of the items is too light or the total weight of the sample is too light, accuracy cannot be guaranteed. You will get an error message, but piece counting will still be allowed. Consult Table 4-1 or Table 4-2 for minimum piece weights and sample weights.

1. If the items you will be counting require a container, you must first tare the container off by pressing the ZERO key.
2. Press the SAMPLE key. The scale will display "5 0 pcs". The scale is prompting you to place five identical items on the platform.

NOTE: If you wish to change the sample number, simply press the SAMPLE key repeatedly until the desired sample number appears. Available choices are 5, 10, 20, 50, 100, 200, and 500.

3. Place the sample items on the platform all at once. Once the scale stabilizes, the scale will display "5 – pcs".
4. Press the SAMPLE key to take the sample. If the sample meets the limits shown in Table 4-1 or Table 4-2, the scale will now display the number of pieces on the scale. If it does not, the scale briefly displays an "Lo", but still allows piece counting. If this occurs, you should use a higher sample amount to achieve better piece count accuracy.

NOTE: If the scale displays an "Add" message, the unit weight of the items you wish to count or total sample weight is too light for your scale to process at all.

MODEL	Capacity / Graduation	Minimum Piece Weight	Minimum Sample Weight
PBS-300	300x 0.01 g	0.008 g	0.25 g
PBS-600	600x 0.02 g	0.016 g	0.5 g
PBS-1500	1500x 0.05 g	0.04 g	1.25 g
PBS-3000	3000x 0.1 g	0.08 g	2.5 g

TABLE 4-1: PBS Piece Counting Sampling Limits

4.2 CLEARING THE PIECE COUNT

1. To clear the piece count, press the SAMPLE key to take a new sample.

4.3 MEMORY ACCUMULATOR

Your scale comes equipped with a handy memory accumulator, which can be used in conjunction with the piece counting feature. As with a hand-held calculator, the memory accumulator can be added to, displayed and cleared at anytime.

NOTE: This feature can be used in any weighing mode. However, the accumulator memory is automatically cleared if the scale is switched to another unit of measure using the UNITS key.

4.3.1 DISPLAYING THE MEMORY ACCUMULATOR

1. Press the MR key. The scale briefly displays the total value stored in the accumulator but only if the Mem annunciator is lit.

NOTE: If a device is connected to the scale's serial port, then that device will receive an automatic printout of the number of entries made and as well as the total value stored in the accumulator. See Appendix B for more detailed information.

4.3.2 ADDING TO THE MEMORY ACCUMULATOR

1. Press the M+ key. The scale shows "Add", lights the Mem annunciator, and then adds the current value shown on the display to the memory accumulator.

NOTE: If a device is connected to the scale's serial port, then that device will receive an automatic printout of that entry. See Appendix B for more detailed information.

4.3.3 CLEARING THE MEMORY ACCUMULATOR

1. Press the MC key. The scale displays "Clr" to indicate that the memory has been cleared and also extinguishes the Mem annunciator.

4.4 PERCENTAGE MODE

This mode is used to indicate the weight percentage of an item you have placed on the scale's platform relative to 100%. This is useful for filling and moisture determination applications.

The scale uses the sampling method to determine the percentage value of 100%. Consult Table 4-3 and Table 4-4 for acceptance criteria.

1. Press the UNITS key until the scale is in percentage mode, indicated by the % sign. The scale should display "100 0 %". If it does not, press the SAMPLE key.
2. If the items you will be counting require a container, you must first tare the container off by pressing the ZERO key.

3. Place the reference item (100%) on the platform. Once the scale stabilizes, the scale will display “100 – %”.
4. Press the SAMPLE key to take the sample. If the sample meets the limits shown in Table 4-3 or Table 4-4, the scale will now display 100%.

NOTE: If the scale displays an “Add” message, the weight of the reference item you wish to sample is too light for your scale to process at all.

MODEL	Capacity / Graduation	Minimum Sample Weight	Best Sample Weight
PBS-300	300x 0.01 g	1 g	100 g
PBS-600	600x 0.02 g	2 g	200 g
PBS-1500	1500x 0.05 g	5 g	500 g
PBS-3000	3000x 0.1 g	10 g	1000 g

TABLE 4-3: PBS Percentage Sampling Limits

4.5 DISPLAY APW MODE

This mode is used to indicate the Average Piece Weight (APW) of the items you are counting in grams.

1. Sample the items you wish to count by following the procedure described in Section 4.1.
2. Press the UNITS key until the APW annunciator is lit. The scale now briefly displays the current APW before reverting back to weighing mode.

CHAPTER 5: SCALE CONFIGURATION

5.1 SCALE CONFIGURATION OVERVIEW

The scale contains two main setup menus: The User (“A”) menu - which configures the serial communication port and enables some user options - and the Factory Setup (“F”) menu. Both menus consist of several menu selections, each with its own sub-menu of choices.

5.2 USER (“A”) MENU

5.2.1 ENTERING THE USER MENU

1. Turn OFF the scale.
2. Press and hold the UNITS key while turning the scale back ON.
3. When the scale shows “- - - - -”, you may release the UNITS key. The scale shows “A 1”.

5.2.2 NAVIGATING IN THE USER MENU

Use the directional keys shown in Figure 5-1 to move around in the User Menu Chart shown in Figure 5-2 on the following page.

1. To move to a new “A” heading, use the M+ (left) or MR (right) key to move right or left in the User Menu Chart.
2. To move to the selection level, press the SAMPLE (down) key once. The current saved selection is shown.
3. To view the available selections for the current “A” heading, use the M+ (left) or MR (right) key to move through the selection field.
4. To save a new selection, press the PRINT (Save) key. To exit without saving, press the ZERO (up) key to return to the current “A” heading.
5. Repeat Steps 2 through 5 until the User Menu is programmed.

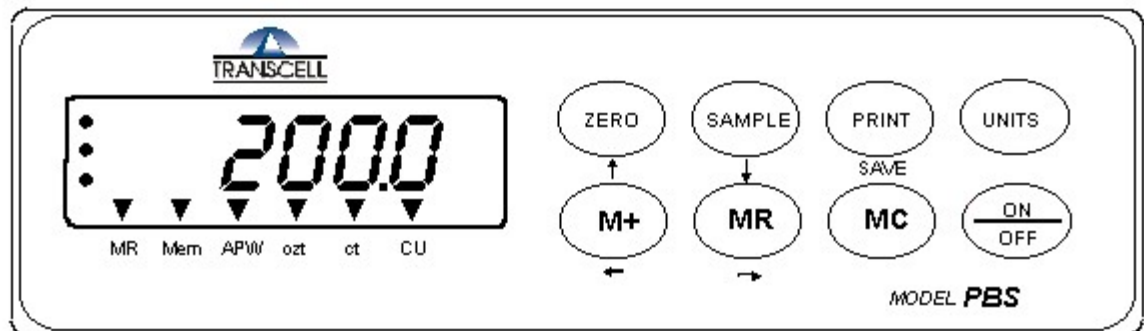


Figure 5-1: User Menu Key Assignments

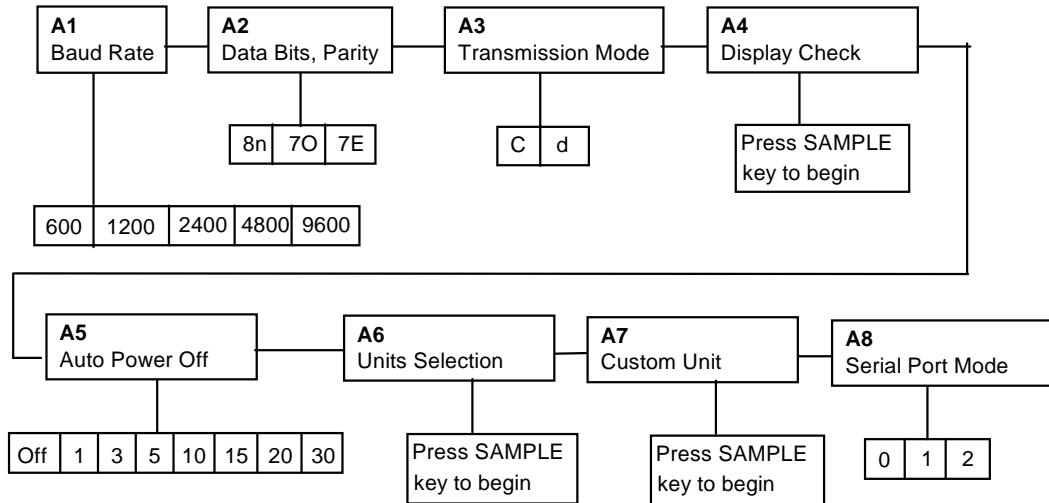


Figure 5-2: User Menu Chart

5.2.3 NOTES ON THE USER MENU

1. Detailed descriptions of the user menu parameters can be found in Table 5-1 on the next page. Factory-set defaults are shown in bold with a checkmark (✓).

NAME/CODE	DESCRIPTION	CODE/VALUE
A1 Baud Rate	Selects the baud rate for data transmission through the serial port.	600 1200 2400 4800 9600 √
A2 Data Bits and Parity	Selects the number of data bits and parity of serial transmission. "8n" = 8 data bits with no parity bit and one stop bit "7O" = 7 data bits with odd parity bit and one stop bit "7E" = 7 data bits with even parity bit and one stop bit	8n √ 7O 7E
A3 Mode of Serial Transmission	Selects when data will be sent out of the serial port to a printer or computer: "C" = Continuous mode; send data continuously "d" = Demand mode; send data when a PRINT command is issued from the computer.	C d √
A4 Display Check	Actuates the function that illuminates all digit segments, decimal points, and LCD annunciators in a test sequence. Pressing the SAMPLE key to scroll down one level begins the test sequence.	Press SAMPLE key to begin sequence
A5 Auto Power Off Period	Selects the auto off time period in minutes. Scale must be idle during this period to shut off automatically. "OFF" = Feature Disabled	OFF √ 1 5 10 15 20 30
A6 Units Selection	Actuates the function that allows customization of the displayed weight units. Pressing the SAMPLE key to scroll down one level begins the sequence.	Press SAMPLE key to begin sequence
A7 Custom Unit	Actuates the function that allows setup of the custom unit. Pressing the SAMPLE key to scroll down one level begins the sequence.	Press SAMPLE key to begin sequence
A8 Serial Port Mode	Selects the mode of the RS-232C serial port: "0" = Full Duplex Mode "1" = APW Transmit Mode	0 √ 1

Table 5-1: User Menu Chart

5.2.4 UNITS SELECTION PROCEDURE (A6)

1. While in the A6.sub-menu, press the SAMPLE key. The scale illuminates the lb annunciator and displays the current status (ON or OFF) of the lb unit.
2. Use the M+ key to toggle the status of the unit from ON to OFF.
3. Use the ZERO key to toggle among the available weight units.
4. Repeat Steps #2 and #3 until you have programmed all units.
5. Press the PRINT key to save all of your new settings and return back to the A6 sub-menu prompt.

5.2.5 CUSTOM UNIT PROCEDURE (A7)

The custom unit multiplier is the conversion factor that you would multiply grams by in order to obtain the custom unit of weight. For example, if you wish to display weight in troy ounces, you would enter in a multiplier value of 0.64301 (1 g = 0.64301 ozt). The scale automatically determines the proper display increment for your custom unit.

1. While in the User Menu, scroll to "A 7", then scroll down once using the SAMPLE key to enter the custom unit multiplier menu.
2. The display will momentarily show "C U" for the custom unit, followed by the current value for the custom unit multiplier with the decimal point.
3. Use the left or right directional keys (shown in Figure 5-1 above) to move the decimal point position to the correct position. Pressing the M+ key or the MR key will change the position of the decimal point.
4. After setting the correct decimal point position, press the PRINT key to save it.
5. The display will momentarily show "SET", followed by a value with one flashing digit. This value will be the current multiplier value with the decimal point position programmed above.
6. Use the four directional keys (shown in Figure 5-1 above) to adjust the displayed value to the actual custom unit multiplier value. Increase the flashing digit by pressing the ZERO key. Decrease the flashing digit by pressing the SAMPLE key. Pressing the M+ key or the MR key will change the position of the flashing digit.
7. After setting the exact value, press the PRINT key to save the custom unit multiplier value. The display will show "EndCU" momentarily, then revert back up to A7.

5.2.6 EXITING THE USER MENU


1. Exit the User ("A") menu by first turning the scale OFF, then turning it back ON without holding down any keys. The display will go through a digit check, then settle into Normal Operating mode. All front panel keys will now return to their normal mode of operation.

5.3 FACTORY SETUP ("F") MENU

To enter the "F" menu, power down the scale. Press and hold the PRINT key while powering on the scale until the scale displays a row of dashes (-----) and beeps. Release the PRINT key. Scale now displays F1. Use the left and right arrow keys to move to another F number. Use the down arrow key to enter the F number. Use the left and right arrow keys to view and select the choices. Press SAVE button to accept choice. Use up arrow key to return to F menu.

To exit F menu, power off scale. Power on scale.

Refer to the "F" menu flow diagram and table for more information.

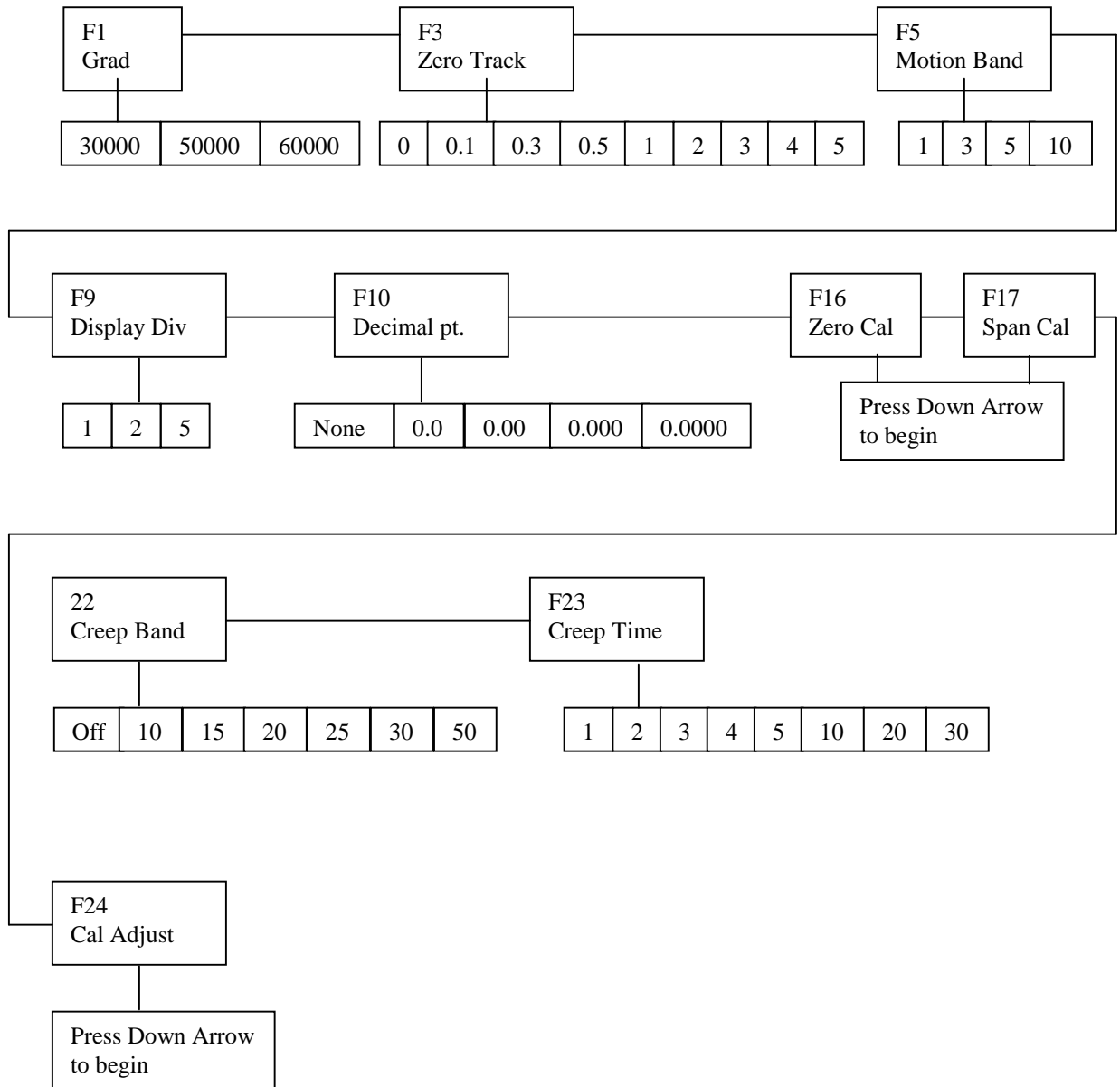


“F” Menu Flow Chart

PARAMETER (Soft Key Text)	DESCRIPTION	AVAILABLE CHOICES
F1 Graduations	Specifies number of full-scale graduations. Value should be consistent with legal requirements and environmental limits on the useful system resolution.	1,000 2,000 5,000 10,000 24,000 30,000 50,000
F3 Zero Track	Selects the range within which the scale will automatically zero. Note that the scale must be in standstill to automatically zero. Selections are in Display Divisions.	0d 0.5d √ 1d 3d 5d
F5 Motion Band	Sets the level at which motion is detected by comparing the present display update with the previous one. If motion is not detected for two seconds or more, scale is in standstill and can process a Print or Zero command. Maximum value varies depending on local regulations.	1d √ 3d 5d 10d
F9 Display Divisions	Determines the desired weight increments. Value should be consistent with legal requirements.	1 2 5
F10 Decimal Pt.	Determines location of the decimal point.	0 0.0 0.00 0.000 0.0000
F16 Zero Calibration	Performs zero calibration. <u>See chapter 6 for calibration instructions</u>	
F17 Span Calibration	Performs scale calibration. <u>See chapter 6 for calibration instructions</u>	
F22 Creep Tracking (CrpTrk)	Sets the range within which the scale will attempt to offset the effects of load cell creep. Use a lower setting for lower capacity load cells and vice-versa. Selections are in seconds. FACTORY SET – DO NOT CHANGE	Off 10, 15, 20, 25, 30, 50,
F23 Creep Time	Sets the range within which the scale will attempt to offset the effects of load cell creep. Use a lower setting for lower capacity load cells and vice-versa. Selections are in seconds. FACTORY SET – DO NOT CHANGE	1, 2, 3, 4, 5, 10, 20, 30,
F24 Calibration Adjust	Allows experienced service technicians to adjust calibration to compensate for linearity adjustments. USE WITH CAUTION	

CHAPTER 6: SCALE CALIBRATION

6.1 SCALE CALIBRATION

1. Enter F menu by powering down scale. Press and hold the PRINT key while powering on the scale until the scale displays a row of dashes (-----) and beeps. Release the PRINT key. Scale now displays F1.
2. Use the RIGHT ARROW key to move to F16 and press the DOWN ARROW key. Scale displays C0 and then displays the zero counts. Press the ZERO key. Scale displays ZERO then 0. Press the SAVE key to accept and return to F16.
3. Use RIGHT ARROW key to go to F17. Use DOWN ARROW key to begin calibration. Scale displays C1 then displays 0.00 with the right digit blinking. Use the LEFT ARROW key to position the blinking cursor over the first digit of the test weight value. Use the UP ARROW key to increment the digit to the test weight value. Place the test weight on the scale. The left most digit blinks faster then slows. After weight has stabilized, press the SAVE key to accept the first cal point.
4. The scale displays SAVE, then SET then C2. Repeat step 3 for cal point 2 and cal point three.
5. To exit without performing three point calibration, press SAVE twice.
6. Scale returns to F17.

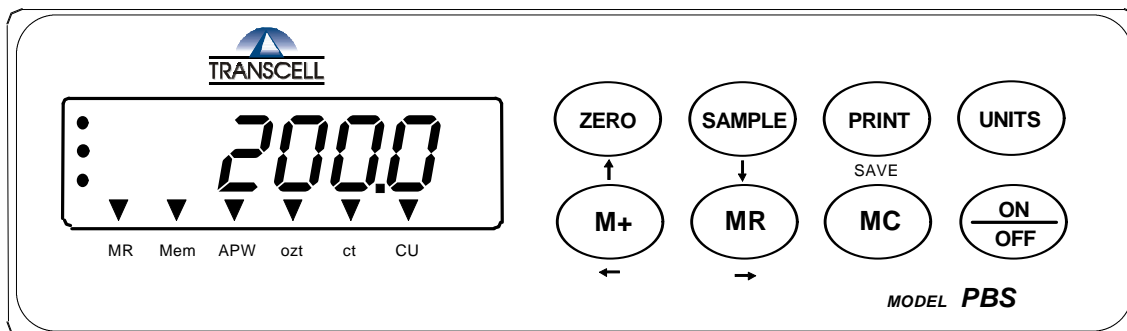

Model	Capacity/Graduation	Test Weight
PBS-300	300 g x 0.01 g	100g
PBS-600	600 g x 0.02 g	100g
PBS-1500	1500 g x 0.05 g	500g
PBS-3000	3000 g x 0.1 g	500g

TABLE 6.1: PBS Calibration Test Weights

6.2 USING CALIBRATION ADJUSTMENT

1. This is an *after scale calibration function*. Be sure to have completed either single or multi point calibration procedures before proceeding. This is not a substitution for scale calibration but is a fine tuning tool to make small compensation adjustments.
2. Experienced service technicians may use the calibration adjustment feature to make fine adjustments for critical applications.
3. Refer the to the “F” menu procedures and use F24.
4. Press the DOWN ARROW key.
5. The scale shows a regular scale display.
6. Place a test weight on the scale and view the results.
7. If the weight is slightly off, press the PRINT key and the right most digit blinks.
8. Use the UP or DOWN arrow key to adjust the digit.
9. Press SAVE to store and return to F24.
10. Repeat procedure with additional weight to adjust linearity.

APPENDIX A: SPECIFICATIONS

CONSTRUCTION:

Housings: Grey ABS
Sub-Platform: Metal
Platter: Plastic with Stainless Steel Cover
Feet: Adjustable Non-skid Hard Rubber

DISPLAY:

6 digit, 7-segment LCD

KEYPAD:

8-key Tactile Keypad

OVER CAPACITY ANNUNCIATION:

103% of Full Scale Capacity

OPERATING TEMPERATURE RANGE:

32°F to 104°F
(0°C to 40°C)

POWER SOURCE:

AC Adapter, 9VDC, 500 mA,
included

ALTERNATIVE POWER SOURCE:

6-AA alkaline batteries, replaceable

SERIAL PORT:


Full Duplex RS-232,
Female DSUB9

WEIGHT:



Net Weight: 3.0 lb (1.4 kg)
Shipping Weight: 4.8 lb (2.2 kg)

PHYSICAL DIMENSIONS:

PBS-300, PBS-600 and PBSP-500

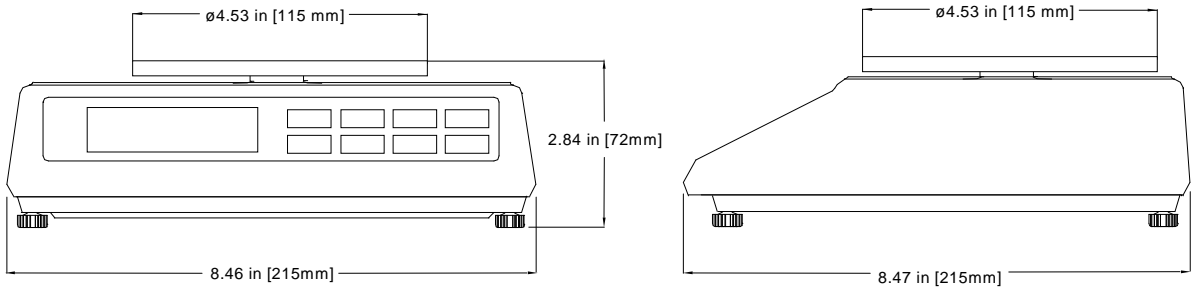

PBS-1500, PBS-3000 and PBSP-5000

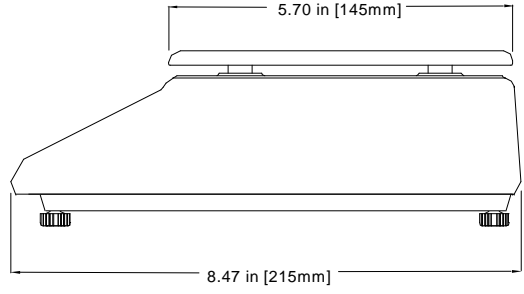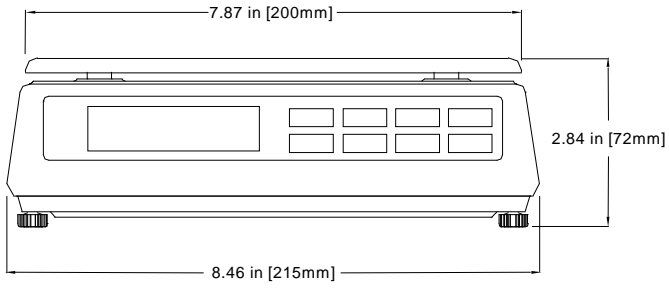

APPENDIX B: SERIAL PORT INFORMATION

B.1 SERIAL PORT HARDWARE

The COM1 serial port is a full duplex RS-232 port designed for connection to one of the following devices:

- Serial receipt printer (dot matrix)
- Computer
- Remote display
- TC-200, TC-2001, TC-2005 or TC-2001 Scale (to transmit APW)

Figure B-1 shows the serial port pin assignments for the scale's female DSUB9 connector. Refer to Figures B-2 and B-3 for some suggested cable diagrams.

1. Plug the serial printer or computer communication cable (not included) directly into the scale's DSUB9 serial port connector.

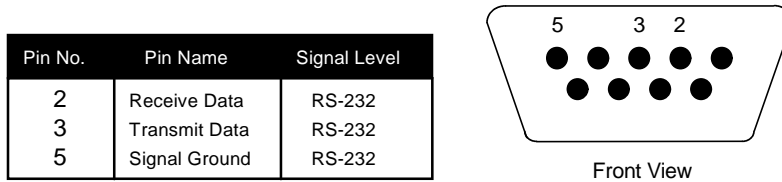


Figure B-1: Pin assignments for the DSUB9 serial port connector

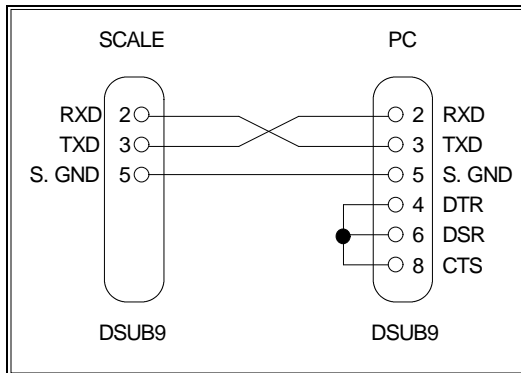


FIGURE B-2. Cable Diagram for Scale to PC

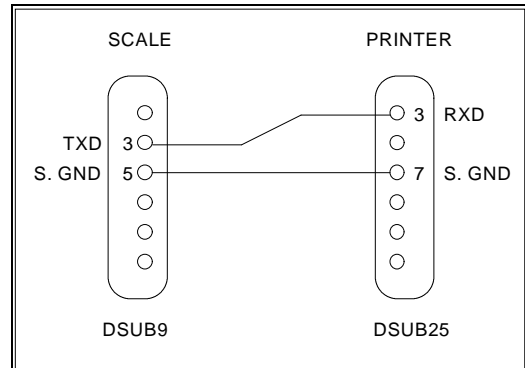


FIGURE B-3. Cable Diagram for Scale to Printer

B.2 SERIAL PORT MODES

B.2.1 DEMAND MODE

The scale provides a full duplex, demand serial transmission mode. This mode is configured in the A3 sub-menu of the User Menu. See Chapter 5 for more information. The demand mode allows control from a host device, such as a PC, and can also be activated by pressing the PRINT key on the scale's front panel. Figure B-2 shows a suggested cable diagram for interface to a PC.

Demand mode is also used when interfacing to a serial printer. Figure B-3 shows a suggested cable diagram for interface to a serial printer.

Figure B-4 shows the serial output data format for demand mode when a PRINT command is issued to the scale.

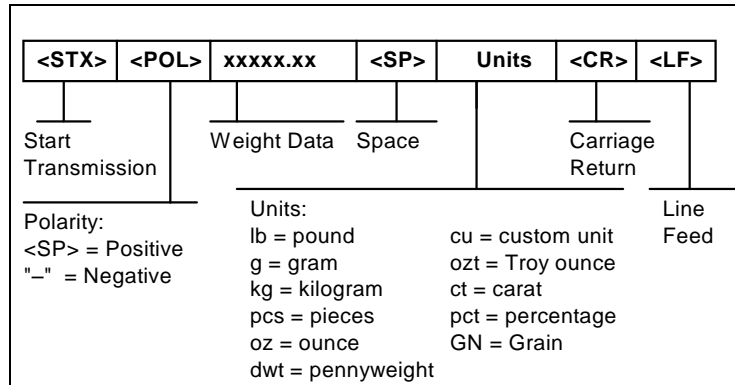


FIGURE B-4. Demand Mode Output Format for Print Command

Figure B-5 shows the serial output data format for demand mode when an M+ (Memory Plus) command is issued to the scale.

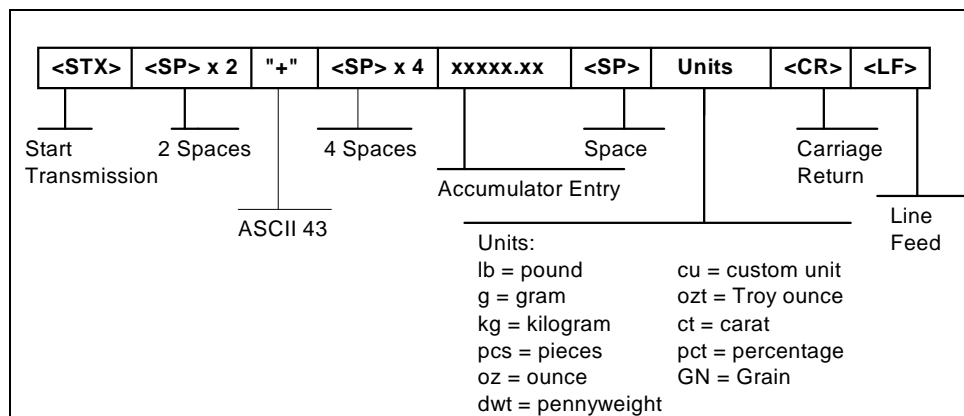


FIGURE B-5. Demand Mode Output Format for M+ Command

Figure B-6 shows the serial output data format for demand mode when an MR (Memory Recall) command is issued to the scale.

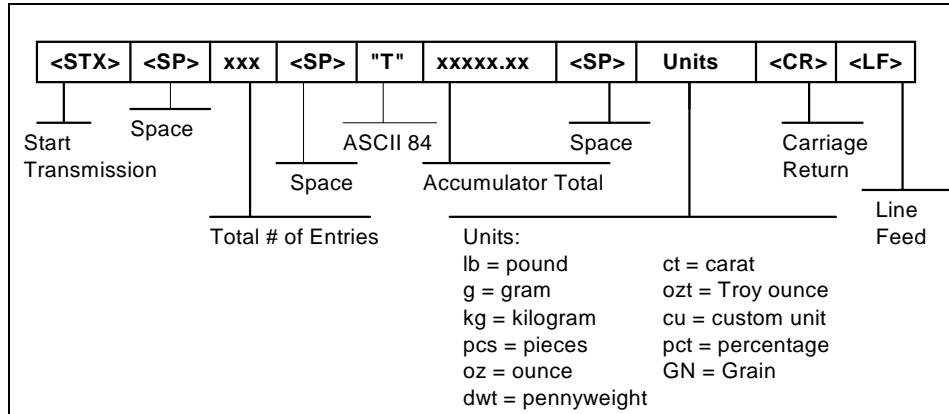


FIGURE B-6. Demand Mode Output Format for MR Command

B.2.2 CONTINUOUS MODE

The scale provides a full duplex, continuous serial transmission mode. This mode is configured in the A3 sub-menu of the User Menu. See Chapter 5 for more information. The continuous mode is used to interface to computers, scoreboards, and other remote devices requiring constant data updating. The transmission occurs at the end of each display update. Figure B-7 shows the serial data format for Continuous Mode.

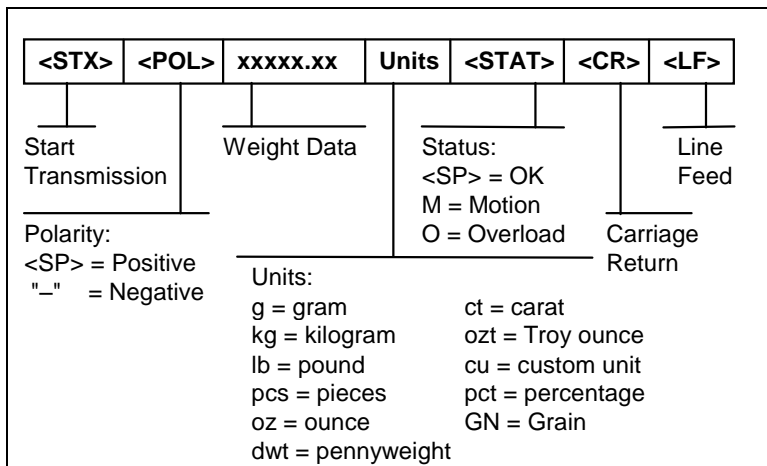


FIGURE B-7. Continuous Mode Output Format

B.2.3 RECOGNIZED HOST COMMANDS

- “P”** - This command is sent to the scale to print the indicated display. The scale will not respond if the scale is in motion, positive overload or negative overload.
- “Z”** - This command is sent to the scale to zero the scale. The scale will not respond if the scale is in motion, positive overload or negative overload.
- “M”** - This command is sent to the scale to add an entry to the accumulator (M+).
- “R”** - This command is sent to the scale to recall the contents of the accumulator (MR).
- “C”** - This command is sent to the scale to clear the contents of the accumulator (MC).
- “U”** - This command is sent to the scale to toggle among the configured units.

B.3 TRANSMIT APW MODE

The Transmit APW Mode is designed specifically for connection to any of the following TC Series scale models:

- TC-200
- TC-2001
- TC-2005
- TC-2010

Its purpose is to allow APW's derived on the PBS precision balance to be transmitted to a larger scale for actual piece counting there. Just press the PRINT key at any time on the PBS to initiate the transmission. Consult your TC-200, TC-2001, TC-2005 or TC-2010 scale manual for more information.

APPENDIX C: ERROR MESSAGES

C.1 ERROR MESSAGES

If the scale encounters an error condition, it will display a message alerting the operator. A description of each display follows:

C.1.1 OPERATOR ERRORS

Message	Explanation
<i>“o o o o o o”</i>	Indicates that the weighing capacity of the scale has been exceeded.
<i>Add</i>	Indicates that there is not enough internal resolution to calculate the unit weight of an item. This means that the items you are counting are too light for the scale to process at all.
<i>Lo</i>	Indicates that unit weight of the items you are sampling is too light for the scale to process accurately.
<i>O L</i>	Indicates that value in the accumulator exceeds the display capacity of the scale.
Err 10	Scale cannot read serial number. Scale must be returned for service.
Err 11	Serial number mismatch. Re-enter setting code from product ID tag.
Err 12	Invalid setting code. Re-enter proper setting code from product ID tag.

Note: Contact Service Department at Transcell for instructions on how to re-enter setting code.